

DNA Stain Type

Ethidium
Bromide

SmartGlow™
SYBR® Green

Excitation Source	DNA Stain Type	
	Ethidium Bromide	SmartGlow™ SYBR® Green
	SmartDoc blue illumination base	Not recommended
Blue light transilluminator	E5001-ORANGE Orange PMMA Filter	E5001-ORANGE Orange PMMA Filter
UV transilluminator	E5001-590 590nm Filter	E5001-535 535nm Filter

For optimum image quality:

- Remove the phone's case
- Line up the camera lens with the photo filter

Place the photo filter into the recess on the top platform

Remove the extender for close ups or small gels



E5001-ORANGE
Orange PMMA Filter
(use with blue light)



E5001-535
535nm Filter
(use with UV)



E5001-590
535nm Filter
(use with UV)